


Instructions for Use

DNAgard® Saliva is designed for the collection of DNA from human saliva samples.

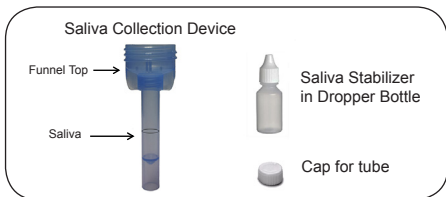

Read all instructions prior to collection.

Precaution

Rinse mouth with water 30 minutes prior to providing a saliva sample, and do NOT eat or drink until finished with saliva collection.

Procedure

1. Remove cap from the dropper bottle and set aside. Spit into the Saliva Collection Device until your saliva (not just the bubbles) has reached the "fill" level marked on the tube. This may take a few minutes.
2. Squeeze the dropper bottle to empty the contents into the Saliva Collection Device.
3. Unscrew the funnel top from the Saliva Collection Device and discard. Be careful not to spill any of the liquid.
4. Screw the white screw cap onto the Saliva Collection Device until it is tightly closed.
5. Mix gently by inverting the Saliva Collection Device 5 to 7 times. The saliva mixture is now ready to ship or store.

Shelf Life

DNAgard Saliva can be stored at room temperature for up to 30 months prior to use.

Product Use Limitations

For Research Use Only. Not for use in diagnostic procedures.

Safety Information

This product is intended solely for the safe collection of saliva samples. Do NOT ingest DNAgard Saliva Solution. If this solution comes in contact with eyes or skin, wash affected areas thoroughly with water. See MSDS at www.biomatrica.com.

Technical Assistance

If you have any questions, please contact Biomatrica technical support at:

Phone: USA (866) DRY-MTRX or (866) 379-6879

Web: www.biomatrica.com

Email: info@biomatrica.com

Rinse mouth with water 30 minutes prior to providing a saliva sample, and do NOT eat or drink until finished with saliva collection.

