


Instructions for Use

SalivaGard™ DNA is designed for the collection and preservation of human saliva DNA.

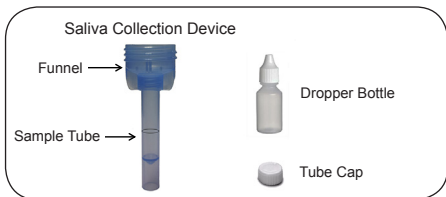

Read all instructions prior to collection.

Precaution

Rinse mouth with water 30 minutes prior to providing a saliva sample. Do NOT eat, drink, smoke, or chew gum until finished with saliva collection.

Procedure

1. Hold Sample Tube upright. Spit into the Funnel until saliva level (excluding bubbles/foam) has reached the "fill" level marked on the Sample Tube. This may take a few minutes.
2. Remove cap from the Dropper Bottle, and empty the contents into the Saliva Collection Device.
3. Unscrew the Funnel from the Sample Tube and discard. Be careful not to spill any of the liquid.
4. Screw the Tube Cap onto the Sample Tube until it is tightly sealed.
5. Mix gently by inverting the Sample Tube 5 to 7 times. The saliva mixture is now ready to ship or store.

Shelf Life

SalivaGard DNA can be stored at room temperature for up to 12 months prior to use.

Product Use Limitations

For Research Use Only. Not for use in diagnostic procedures.

Safety Information

This product is intended solely for the safe collection of saliva samples. Do NOT ingest SalivaGard DNA solution. If this solution comes in contact with eyes or skin, wash affected areas thoroughly with water. See SDS at www.biomatrica.com.

Technical Assistance

If you have any questions, please contact Biomatrica technical support at:

Phone: 858-550-0308 or 866-379-6879

Email: info@biomatrica.com

Web: www.biomatrica.com

Rinse mouth with water 30 minutes prior to providing a saliva sample, and do NOT eat or drink until finished with saliva collection.

